SOP: Isofluranstation

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SOP: Isoflurane station

General information

The isoflurane station is used for the narcotization of mice in the context of officially approved interventions on laboratory rodents.

Technical settings

- O₂-flow: 0,5 l/min
- Isoflurane concentration: 4 % for initiation, 1 - 2 % for maintenance

Procedure

- Check connections of Isoflurane station
- Check isoflurane level, top up if necessary
- Check the weight of the carbon filters, change filter if necessary (max. + 50g)

a) Anesthesia induction / short interventions:

- Set switch to „Box“
- Turn on O₂ wall tap
- Turn on O₂ flow on isoflurane station
- Set isoflurane flow to initiation concentration
- Wait until the anesthesia box has been filled with isoflurane
- Place mouse in anesthesia box and wait for unconsciousness

b) Longer interventions with anesthesia mask:

- Set switch to „Mask“ and reduce isoflurane concentration to maintenance dosage
- Transfer of mouse to anesthesia mask

After completion of the intervention: Switch off the isoflurane and O₂ flows; turn off the O₂ wall tap

Measures for spillage of isoflurane (see also operating instructions for isoflurane)

The room must be vacated immediately. An isoflurane emergency kit (mask and chemical binder) is available to remove the spilled isoflurane. Entering the room is only permitted with the full protective mask, spilled isoflurane is absorbed with the help of the chemical binder (granulate). After use, this is placed under the chemical fume hood until the isoflurane has completely evaporated (> 24h). The granulate is then disposed of in the waste for chemically contaminated materials.